



Separation of DTA, TTA and CTA

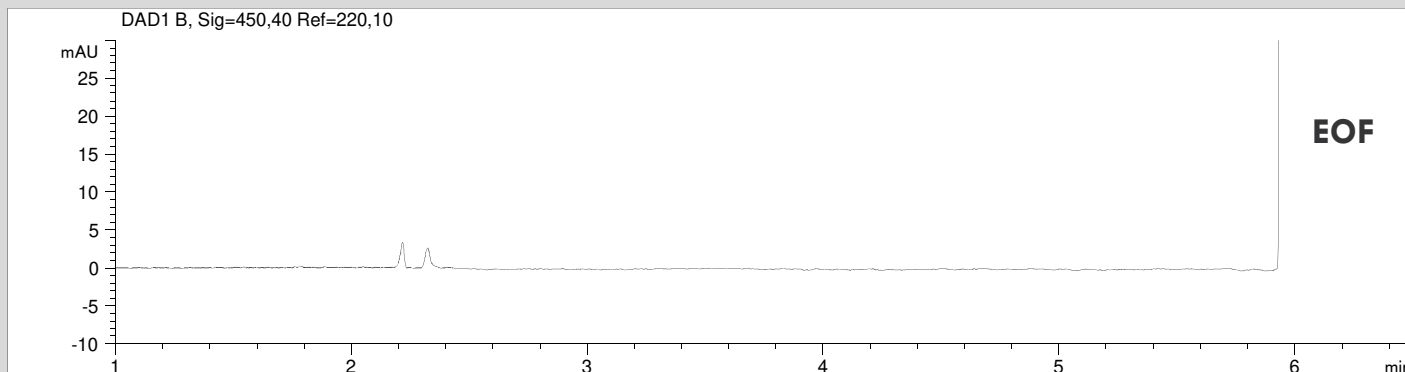

Blank solution: pure water

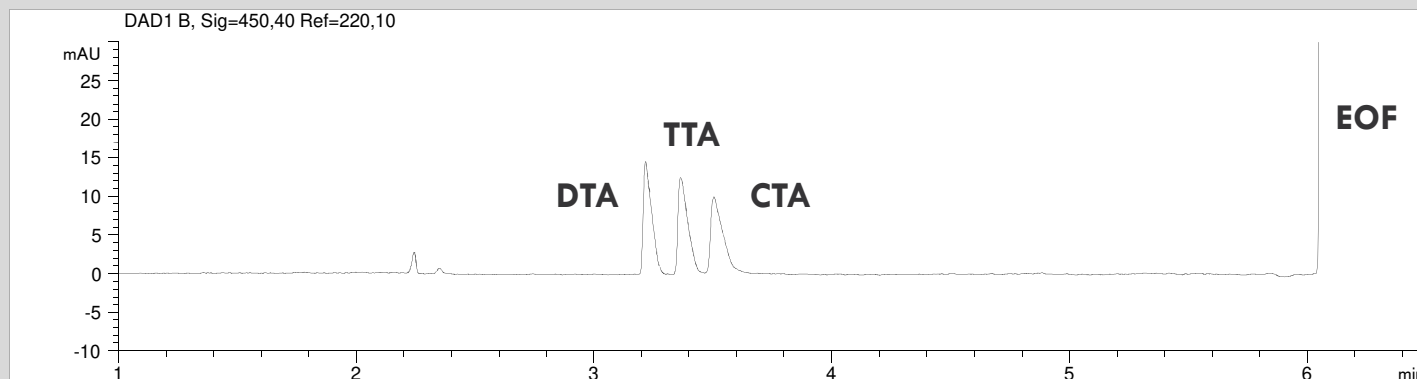

Standard mix solution: Dodecyltrimethylammonium cation, Tetradecyltrimethylammonium cation and Cetyltrimethylammonium cation, 50 mg/l each

- **Mode: CZE**
- **Electrolyte: ABI**
- **Capillary: 50 μ m ID, 48 cm total**
- **Separation: +30 kV, 25°C**
- **Injection: hydrodynamic, 50 mbar, 20 s**
- **Detection: indirect UV, 220 nm**