




## Determination of vitamin C in food supplements

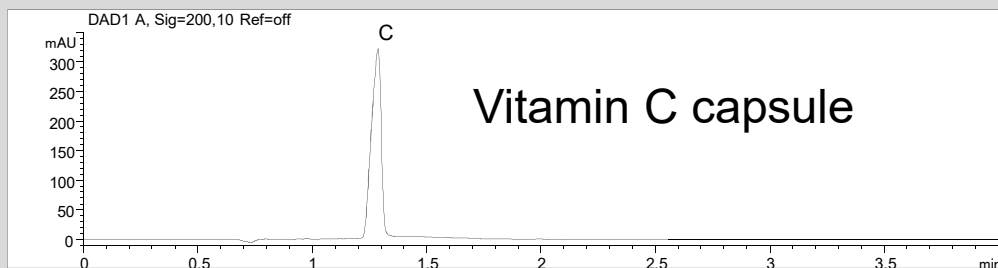

Sample preparation: 100 mg / 10 ml 1:5

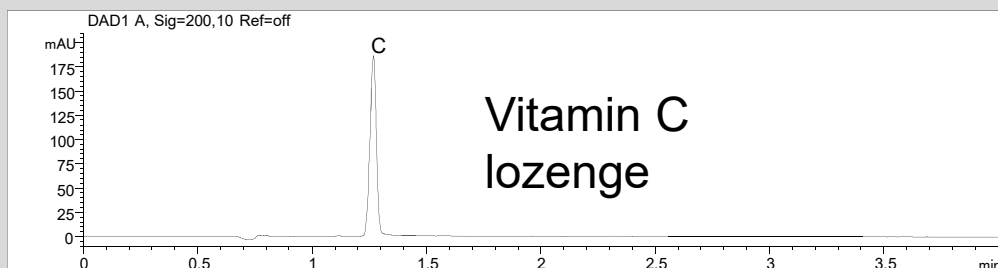

Sample preparation: One tablet in 20 ml, dilution 1:10

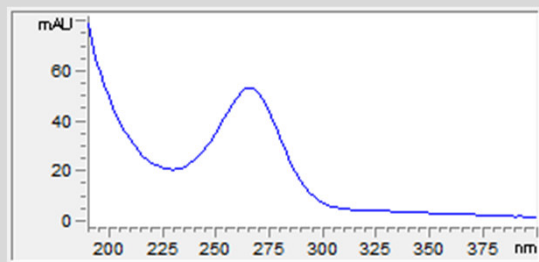

- **Mode: MEKC**
- **Electrolyte: Borate/SDS**
- **Capillary: bubble cell, 50 µm ID, 56 cm**
- **Separation: 30 kV, 25°C**
- **Injection: hydrodynamic, 50 mbar, 20 s**
- **Detection: direct UV, 200 nm**

Operator: C. Amoussou, Chr. Öffler, 07-2021