



Determination of anions and organic acids in food supplements

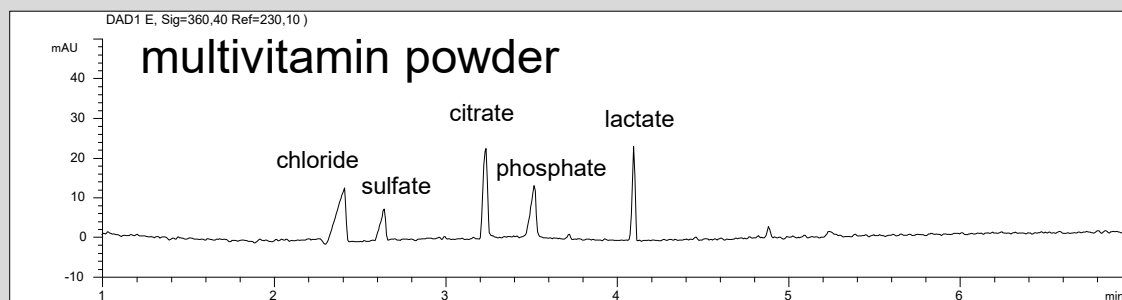

Sample with chloride, sulfate, citrate, phosphate, lactate

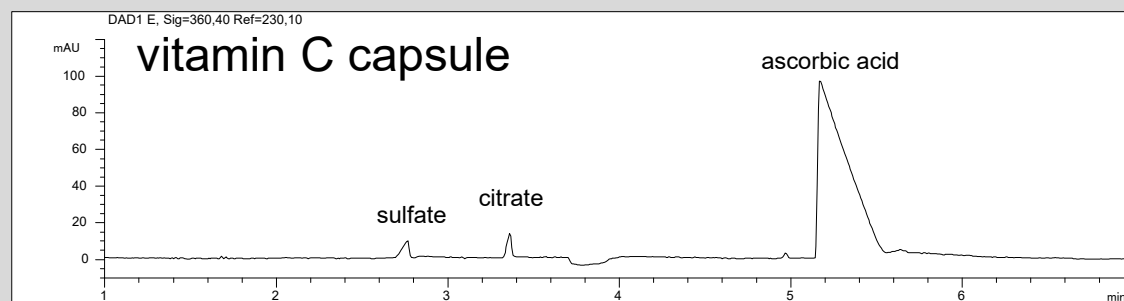

Sample with sulfate, citrate, ascorbic acid

- **Mode: CZE**
- **Electrolyte: salicylate**
- **Capillary: 50 µm ID, 64 cm total**
- **Separation: -30 kV, 25°C**
- **Injection: hydrodynamic, 50 mbar, 10 s**
- **Detection: indirect UV, 230 nm**

Operator: C. Amoussou, Chr. Öffler, 07-2021